

Specification Sheet

CELLINK Bioink

Product description	CELLINK Bioink is the first universal bioink designed to 3D bioprint human tissue models. It is a hydrogel which exhibits an excellent printability and is composed of alginate and highly hydrated cellulose nanofibrils which supports 3D cell culture of various cell types. CELLINK Bioink crosslinks easily with the included Crosslinking Agent. Culture in medium supplemented with calcium to maintain the structure for longer culture periods. For description on how to mix with cells, bioprint and crosslink, follow the Bioprinting Protocol .
Intended use	Biocompatible material for 3D bioprinting, <i>Research Grade</i> . For research use ONLY. Not intended for <i>in vitro</i> diagnostics or <i>in vivo</i> uses. Not intended for administration in humans or animals. Produced under sterile and aseptic conditions.
Product number	IKC20000.
Shelf life	Minimum 6 months, expiration date stated on package.
Storage and handling	Store at 4-8°C. DO NOT FREEZE. Avoid temperature fluctuations. Ensure that the bioink container is capped prior to storage to prevent drying.
Safety	Handle in accordance with good hygiene and laboratory safety practices. Read Safety Data Sheet for more information regarding ingredients and potential hazardous compounds.
Related documents	Bioprinting Protocol as well as Safety Data Sheet can be downloaded from our website https://cellink.com/product/cellink-bioink . Scan the QR code below to reach the product webpage



Property of final bioink	Specification	Method
<i>Appearance</i>	White translucent gel	Visual inspection.
<i>Sterility</i>	Sterile	Tested for the presence of bacteria, fungi and yeast. Tested on raw materials.
<i>Endotoxin level</i>	<5 EU/mL	Limulus Amoebocyte Lysate assay, Pharmacopoeia 2.6.14 "Bacterial endotoxins": Method D, accredited by SWEDAC. Accreditation Certification 1240: ISO 15189, 2010-11-22. Tested on raw material components.
<i>Viscosity</i>	$\geq 7 \text{ kPa}\cdot\text{s}$ at 0.01 s^{-1} ; $\leq 3 \text{ Pa}\cdot\text{s}$ at 200 s^{-1}	Tested using rotational 20 mm plate-plate HR-2 TA Instruments Rheometer. Flow sweep parameters: 25°C, shear rate from 0.001 s^{-1} to 500 s^{-1} .
<i>pH</i>	6.5-7.4	Assessed with pH paper.